

Rely+On™ Virkon™


Broad Spectrum
Virucidal Disinfectant

How to prepare and use a Rely+On™ Virkon™ disinfectant solution

Preparing a 500ml 1:100 disinfectant solution with Rely+On™ Virkon™ 5g tablets


STEP 1

Wear appropriate personal protective equipment before preparation and use of disinfectant solutions

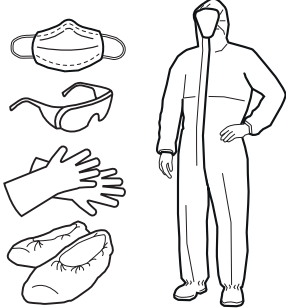

STEP 2

Fill a 500ml trigger spray bottle with room temperature water

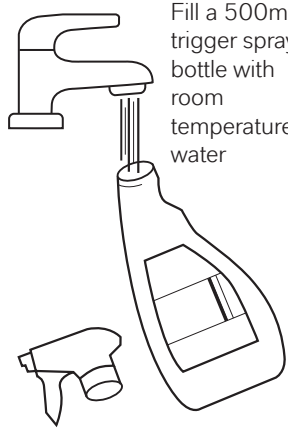


STEP 3

Add 1 x 5g Rely+On™ Virkon™ tablet and wait for it to dissolve completely



STEP 4


The Rely+On™ Virkon™ disinfectant is now ready for use

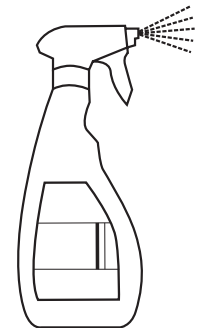

Preparing a 500ml 1:100 disinfectant solution with Rely+On™ Virkon™ powder concentrate


STEP 1

Wear appropriate personal protective equipment before preparation and use of disinfectant solutions

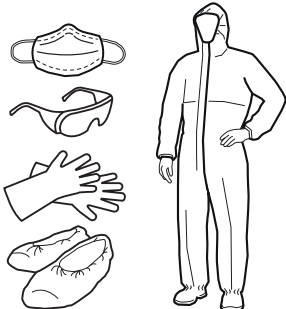


STEP 2

Fill a 500ml trigger spray bottle with room temperature water

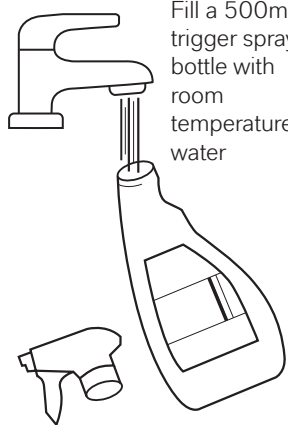

STEP 3

Weigh out 5g Rely+On™ Virkon™ powder concentrate

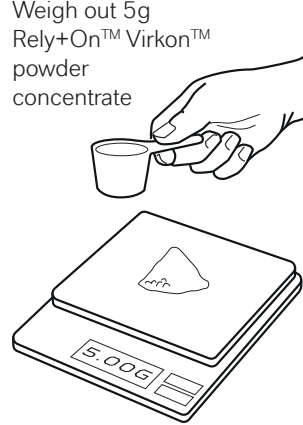

STEP 4

Add the 5g to the bottle and wait for it to dissolve completely



Applying Rely+On™ Virkon™ disinfectant solution via a trigger spray bottle


STEP 1

Spray Rely+On™ Virkon™ solution onto the surface to be disinfected



STEP 2

Wait a minimum of 10 minutes

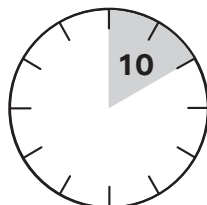


STEP 3

Spray all surfaces that have had disinfectant applied with fresh clean water



STEP 4


Then wipe surface dry with a clean cloth

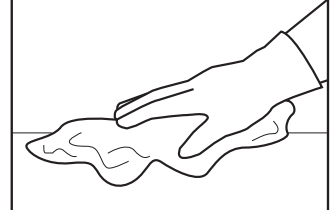

Preparing 1:100 disinfectant solutions with Rely+On™ Virkon™ powder concentrate


STEP 1

Wear appropriate personal protective equipment before preparation and use of disinfectant solutions

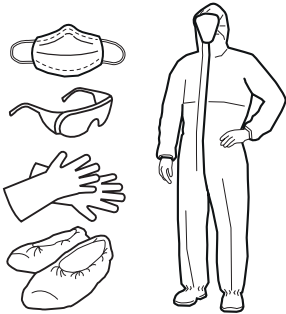


STEP 2

Fill a suitable container with the required volume (e.g. 5 litres) of room temperature water

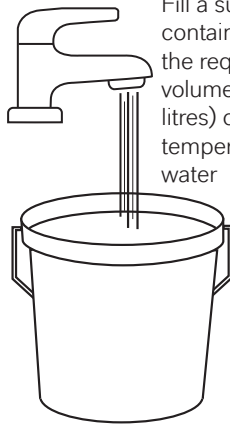


STEP 3

Weigh out 10g of Rely+On™ Virkon™ powder concentrate for every 1 litre of water

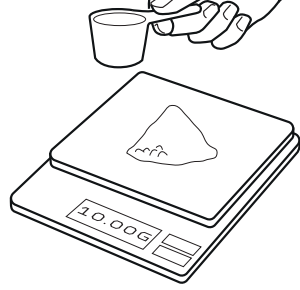

STEP 4

Add the Rely+On™ Virkon™ powder to the water and then stir until completely dissolved. The disinfectant solution is now ready to use

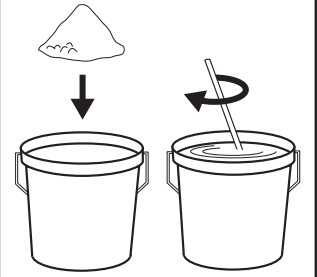

Applying Rely+On™ Virkon™ disinfectant solutions


STEP 1

To disinfect large areas of floor space, apply Rely+On™ Virkon™ solution using a mop and bucket




STEP 2

To disinfect walls, doors, and other large surfaces, use a bucket and sponge or cloth to apply Rely+On™ Virkon™ disinfectant solution

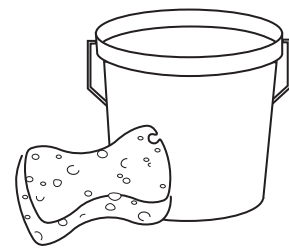

STEP 3

Wait for a minimum of 10 minutes

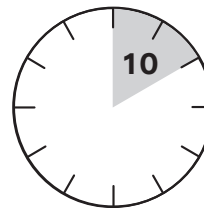

STEP 4

Rinse all surfaces with clean water when materials compatibility is of concern



* Not for use for the disinfection of medical devices.

** Users should always check for compatibility with any sensitive materials before use. Do not use on any soft metals such as brass or copper. Not recommended for use on soft furnishings, textiles or carpets.

The specified uses and registered claims may vary from country to country. Please contact LANXESS to verify country-specific approved uses.

Use biocides safely. Always read the label and product information before use.

LANXESS
Energizing Chemistry

Antec International Limited
Windham Road, Chilton Industrial Estate
Sudbury, Suffolk CO10 2XD,
United Kingdom

Tel: +44(0)1787 377305
relyondisinfection@lanxess.com
www.relyondisinfection.com

©2020 LANXESS. Rely+On™ Virkon™, LANXESS™ and the LANXESS Logo are trademarks of LANXESS Deutschland GmbH or its affiliates. All trademarks are registered in many countries worldwide.